



## planetary ball mill basic information introduction

### Introduction:

Before operating the Laboratory planetary ball mill, please read the following rules carefully.

### 1.Safety

The small Laboratory ball mill operation and maintenance must comply with instructions.

Misusing machine will cause damage to the ball mill and make it stop working.

### 2.Pay attention to using the power:

Before connecting the power, ensure that the voltage is consistent with the machine required voltage (allowing a deviation of +/- 10%), and ensure the rated load of powder socket should be no less than machine requirement.

### 3.Pay attention to using the power cord:

The machine usually uses incidental powder cord. If the power cord is broken, it must be replaced not be repaired. And the same type and specification of the power cord should be replaced. When it is working, it is not allowed to put any items on the power cord. Especially do not put it in the area where people often walk around.

### 4.Pay attention to the machine location:

Cool, ventilated, dry, dust-proof environment

